UVisco

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UV-1600/1800 - V1600/1800

CE

LC-Instru

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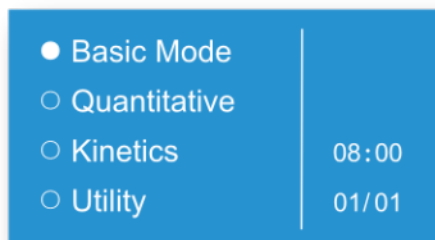
Menu Principal

Move the cursor on the function menu you want, then press ENTER key to go into the corresponding interface.

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Absorbance, Transmittance or Concentration measurements. It can display and save 200 groups of data, 5 groups per screen.

No.	WL.	Abs.
1	230.0	0.001
2	340.0	0.000
3	450.0	0.002
4	540.0	0.000
5	620.0	0.003

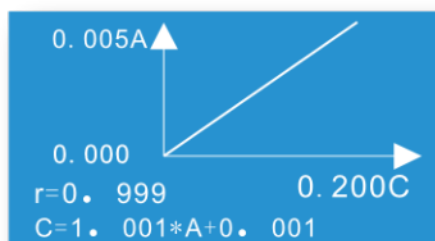
Analyse Quantitative

2.1 Standard Curve

At most 9 standard samples can be used to establish a standard curve. The curve and its equation will be displayed on the screen simultaneously. You can measure your unknown concentration solutions by the curve. Total 100 curves can be saved in the RAM.


2.2 Coefficient Method

If you have known the coefficient k & b of the formula $C=kA+b$, you can input the value directly by the button, then to test your unknown solutions.

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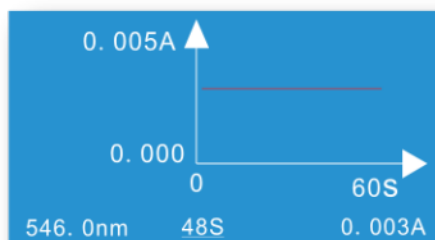This mode may be used for time course scanning or reaction rate calculations. Abs. Vs time graphs is displayed on the screen in real time.

Utilitaires

Lamp management, time & date set, get dark current, calibrate wavelength, default system, and some other system functions.

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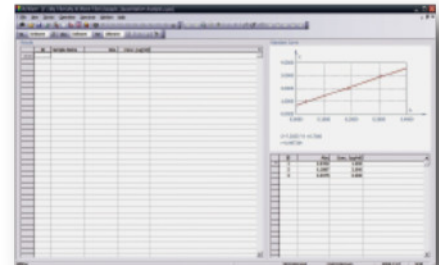
M.Wave Professional application software is based Microsoft Windows®, the instrument can be controlled by PC software through the built-in USB communication port, which makes the UV/V-1 Series with more functions and easy to control.

Analyse Quantitative


Use up to 20 standards to establish standard curve. Three methods for fitting a curve:

- 1-Linear fit
- 2-Linear through zero
- 3-Square fit

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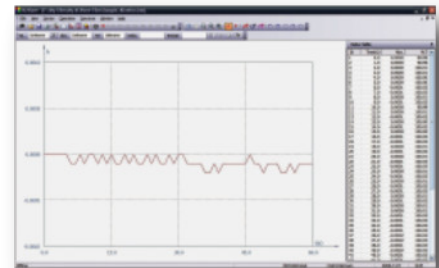The Kinetics mode may be used for time course scanning or reaction rate calculations. Abs. Vs. Time graphs is displayed.


Prise de spectres

Automatically records peaks and valleys. The quantity of the curves stored is unlimited.


Post-run manipulation and processing includes.

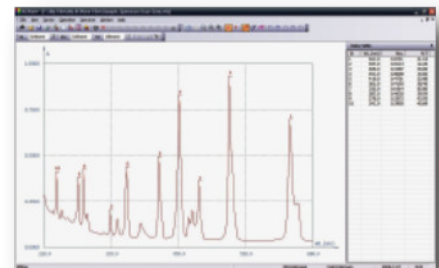1. Re-scaling axes, curve
2. Smoothing, combination, zooming, overlap...
3. 1st to 4th derivative

Analyses multi-longueurs d'ondes

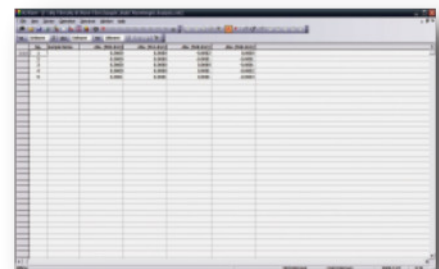You can set up to 20 wavelengths to measure a sample.


Tests ADN/Protéines

Optional two formulas:

$$\text{DNA Concentration} = 62.9 \cdot A_{260} - 36.0 \cdot A_{280}$$

$$\text{Or } 49.1 \cdot A_{260} - 75.8 \cdot A_{230}$$

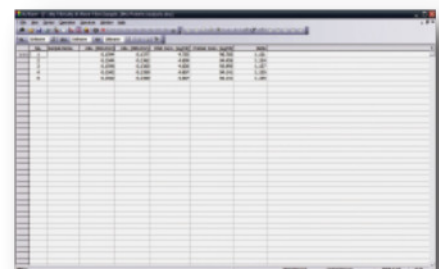You can also enter other wavelengths and factors to calculate.

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